

Animal Control DNAs for LCD-Array kits

Animal Control DNAs are delivered in 0.5 x TE Buffer at 100 ng/μl (+/- 2) and are shipped without cooling (express delivery). Each vial contains 25 μl (2.5 μg in total). Vials should be stored at -20°C upon arrival.

For routine use as positive controls we recommend to dilute the stock solution 1:10 by adding 225 μl 0.5 x TE buffer at arrival and storage of diluted aliquots with 10 ng/μl at -20°C. For sensitivity tests with end dilution rows it is recommended to start from the concentrated stock solution.

Name	Species	Order Code:
Cattle	<i>Bos taurus</i>	ACD-001-025
Sheep	<i>Ovis aries</i>	ACD-003-025
Horse	<i>Equus caballus</i>	ACD-004-025
Goat	<i>Capra hircus</i>	ACD-006-025
Camel	<i>Camelus dromedarius</i>	ACD-007-025
Water Buffalo	<i>Bubalus bubalis</i>	ACD-008-025
Pig, domestic	<i>Sus scrofa</i>	ACD-009-025
Kangaroo	<i>Macropus rufus</i>	ACD-010-025
Hare	<i>Lepus europaeus</i>	ACD-011-025
Rabbit	<i>Oryctolagus cuniculus</i>	ACD-012-025
Reindeer	<i>Rangifer tarandus</i>	ACD-013-025
Roe Deer	<i>Capreolus capreolus</i>	ACD-014-025
Red Deer	<i>Cervus elaphus</i>	ACD-015-025
Fallow Deer	<i>Dama dama</i>	ACD-016-025
Springbok	<i>Antidorcas marsupialis</i>	ACD-017-025
Dog	<i>Canis lupus familiaris</i>	ACD-018-025
Cat	<i>Felis silvestris catus</i>	ACD-019-025
Chicken	<i>Gallus gallus</i>	ACD-020-025
Turkey	<i>Meleagris gallopavo</i>	ACD-021-025
Goose	<i>Anser anser</i>	ACD-022-025
Ostrich	<i>Struthio camelus</i>	ACD-023-025
Mallard Duck	<i>Anas platyrhynchos</i>	ACD-024-025
Muscovy Duck	<i>Cairina moschata</i>	ACD-025-025
European Elk	<i>Alces alces</i>	ACD-027-025

Other species available on request.